

InSpec™

# Technical Data of InSpec Products by Chemistry

## Aqueous Cleaning Detergents



## InSpec™ N7

### Low Toxicology, pH Neutral Detergent Solution

#### Use Instructions

InSpec™ N7 is designed for cleaning hard surfaces by spraying, wiping and mopping. Hold spray approximately 15cm to 20cm from the surface area. 5L versions, pour into an appropriate container for mopping. 100ml concentrate added to 2.5L of water is required for mopping. Apply to surface to ensure complete coverage.

#### Technical Data

N7 Ready to Use (RTU) is a 4% neutral detergent, pH7. The neutral detergent has surfactant and sequestering agents designed to assist the removal of process residue and light soiling in a cleanroom environment.

#### Storage & Safe Handling

Store upright in original closed containers, away from sunlight and extremes of temperature. Full guidance on the handling and disposal of this product is available in the Safety Data Sheet (SDS).

#### Material Compatibility

Application of solutions, when used as directed, will not affect materials normally encountered. Compatibility information in the Technical Dossier.

#### Benefits

- Designed for cleaning floors, walls, ceiling and other hard surfaces for the removal of process derived residue and light soiling.
- Ideal for critical surfaces with soiling near to product manufacturing
- Very low foaming, low residue
- detergent with an excellent low toxicology profile
- Suitable for manual cleaning and available in convenient ready to use and concentrate formats.
- Low concentration
- 3 months validated 'in-use' shelf life
- All formats manufactured with WFI quality water
- Multiple bags for cleanroom transfer
- Manufactured in accordance with GMP

#### Technical User Information

InSpec™ N7 is available in sterile RTU pre-diluted formats and as a concentrate.

InSpec™ N7 is manufactured to GMP in an ISO 6 cleanroom. The product is filled and capped in an ISO class 5 environment.

RTU solutions are diluted with WFI quality water.

The solution is filtered through a 0.2 micron filter and gamma irradiated to a validated process to give a sterility assurance level SAL of 10<sup>-6</sup>

InSpec™ N7 is supplied with Certificates of Conformity, Analysis, Irradiation and Sterility.

For screw cap bottles use the entire contents after opening.

InSpec™ N7 Sterile Products		
Product	Sterile Part No.	No. per case
InSpec N7 1L PREMIUM Trigger Spray	N7WFI1CC-1LS	6
InSpec N7 5L	N7WFI1-5LS	2
InSpec N7 Concentrate - 100ml	N7WFI1-100MLS	50
InSpec N7 Pre-impregnated Wipes x 15	N7WFI1PCH15S	25
InSpec N7 Pre-impregnated Wipes x 50	N7WFI1PCH50S	16

InSpec™ N7 Non-Sterile Products	
Non-Sterile Part No.	No. per case
NSN7WFI-1L	6
NSN7WFI-5L	2
NSN7CONC-100ML	50
N/A	N/A
N/A	N/A

Wipe Media Type	Individual Wipe size (mm)
N/A	N/A
N/A	N/A
N/A	N/A
N/A	N/A
N/A	N/A